



食用油分析

食用榛子油中 10 种甘油三酯的分析

色谱柱: MEGA-LAP 25m x 0.32mm 0.1um df

订货号: C-LAP-032-010-25

MEGA[®]
CAPILLARY COLUMNS
LABORATORY

Triglycerides - Hazelnut Oil
on MEGA-LAP Column

Column: MEGA-LAP - 0.32mm, 0.10 μ m, 25m
Catalog Code: C-LAP-032-010-25

MEGA-LAP (Lipids Analysis Phase) is a dedicated stationary phase optimized for Lipids, Sterols and saturated and unsaturated Triglycerides separations, stable over 370°C.

Conditions:
Injection: Split 300°C, 1.0 μ L, 1:80 Split Ratio.
Detector: FID 300°C.
Oven Temp.: 120°C, 5°C/min, 330°C.
Carrier Gas: Hydrogen, 2mL/min.

Peaks:
1. PPO
2. POS
3. POO
4. PLO + OOPo
5. PLL + PoOl
6. BOO
7. OOO
8. OOL
9. OOL
10. LLL

Acknowledgement: Dott. C. Mariani - Stazione Sperimentale Oli e Grassi, Via Giuseppe Colombo 79 - Milano - Italy.

since 1980
MEGA s.n.c. - Capillary Columns Laboratory - Via Plinio, 29 - 20025 Legnano (MI) - Italy - Tel./Fax: (+39) 0331.54.79.24 - info@mega.mi.it

www.mega.mi.it

